


Quick Reference Card

For safety and biohazard guidelines, refer to the “Safety” section in the *PrepMan® Ultra Sample Preparation Reagent Protocol (PN 4367554)*. For all chemicals in **bold** type below, read the MSDS and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.


For a list of required and recommended equipment, consumables, and reagents refer to the *PrepMan® Ultra Sample Preparation Reagent Protocol (PN 4367554)*.

DNA Sample Preparation for Bacterial and Fungal Cultures

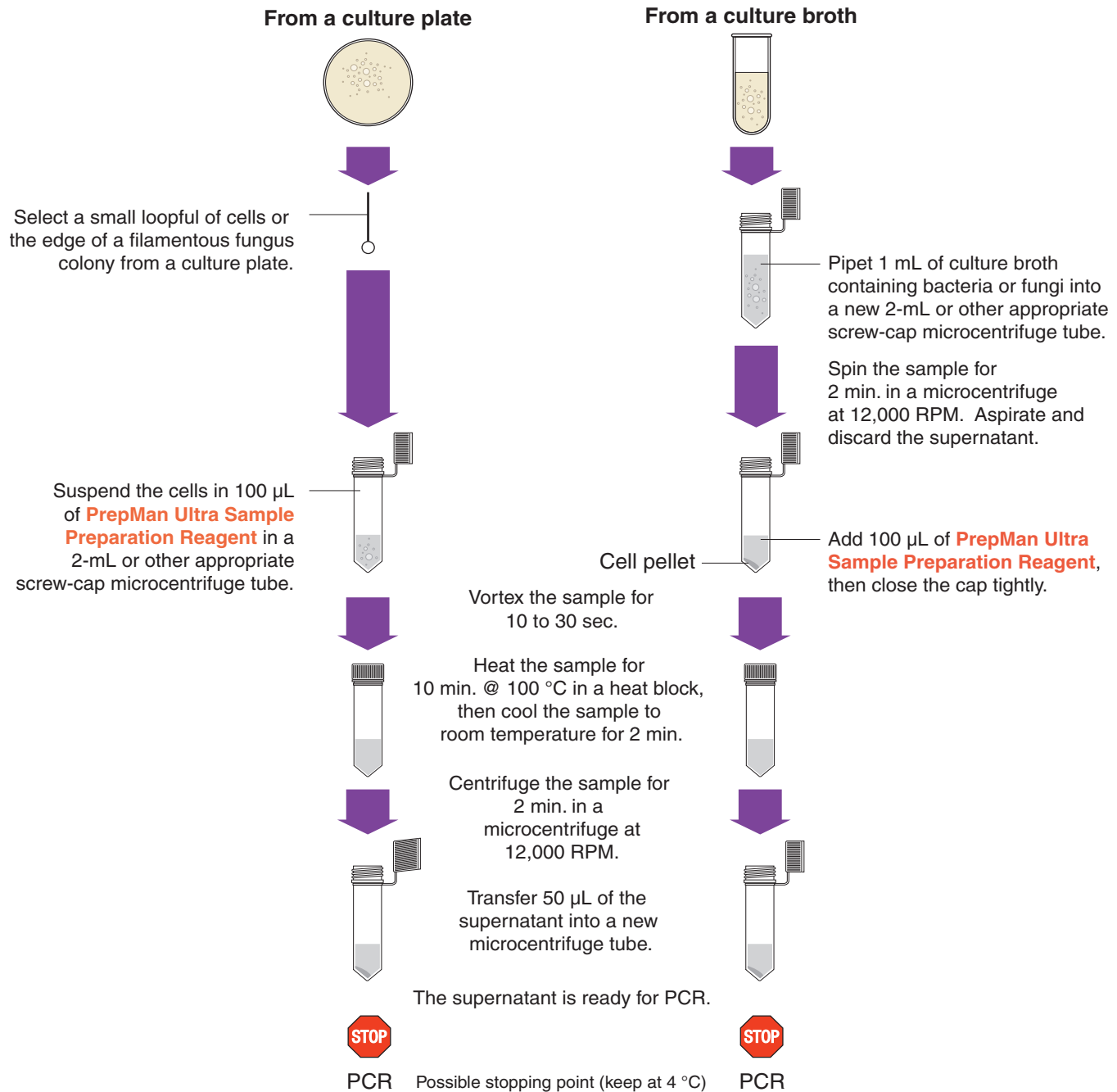

DNA Sample Preparation for Processed Food Samples (GMO detection)

From sample or concentration reference standard

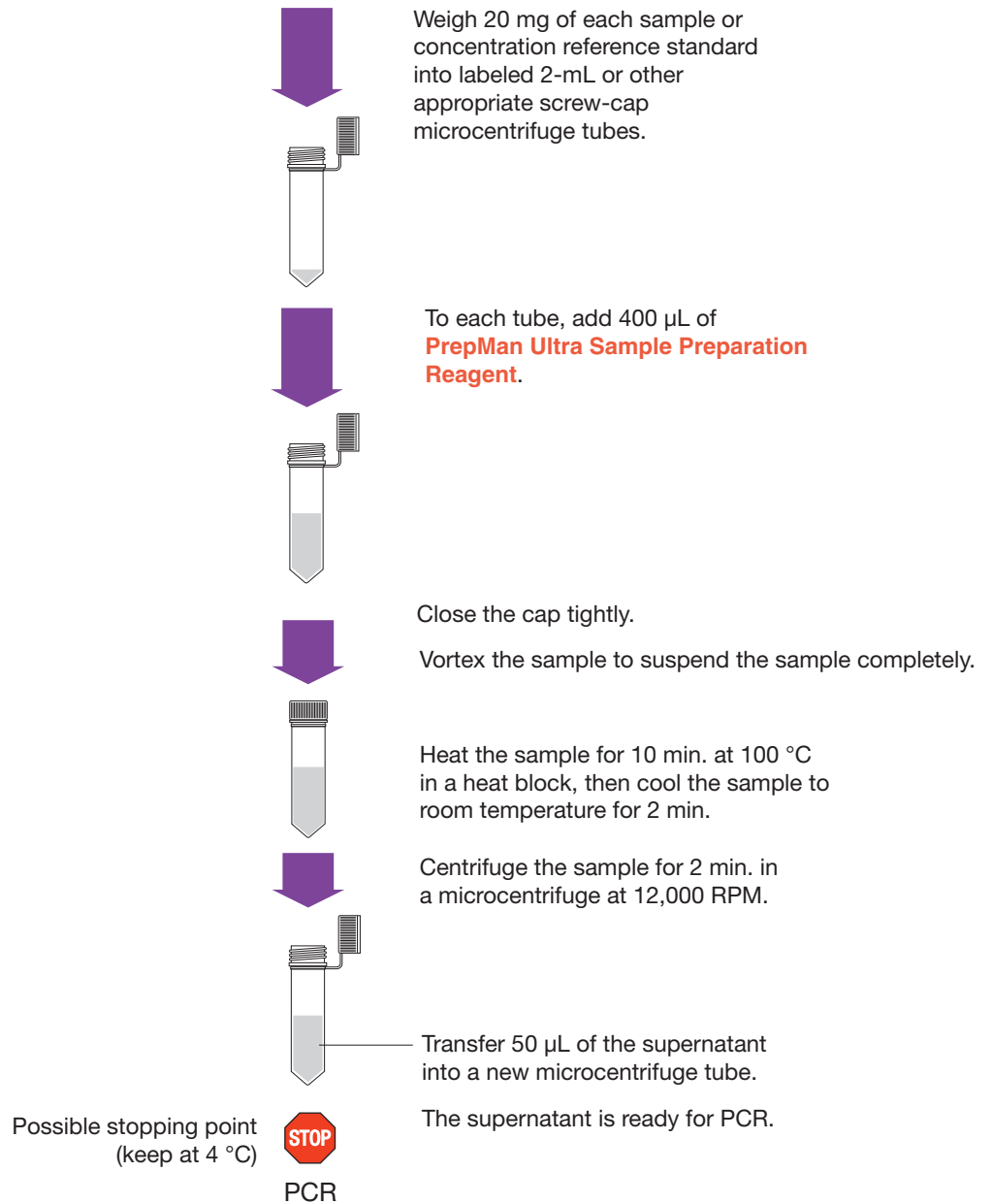

Copyright© 2008. Applied Biosystems. All rights reserved.

For Research Use Only. Not for use in diagnostic procedures.

NOTICE TO PURCHASER:

PLEASE REFER TO THE PREPMan ULTRA SAMPLE PREPARATION REAGENT PROTOCOL FOR LIMITED LABEL LICENSE OR DISCLAIMER INFORMATION.

Applied Biosystems, AB (Design) and PrepMan are registered trademarks of Applied Biosystems or its subsidiaries in the US and or certain other countries.

All other trademarks are the sole property of their respective owners.

03/2008

 Applied Biosystems